



# DURACLONE RE

THERE ARE EVENTS YOU CANNOT AFFORD TO MISS



Your clinical research trial companion


**VISUALIZE**  
*the possibilities.*

For Research Use Only - Not for use in Diagnostic procedures

 **BECKMAN  
COULTER**  
*Life Sciences*

# THERE ARE EVENTS YOU CANNOT AFFORD TO MISS

Rare event cytometry has emerged as a powerful and important tool in clinical research of blood disorders. Sensitive and specific detection of just a few dozens of abnormal target cells among millions of normal cells is one of the most challenging applications in flow cytometry. The DuraClone RE kits have been optimized to accommodate either large sample volume or high cellular concentrations to achieve high sensitivity for rare populations.

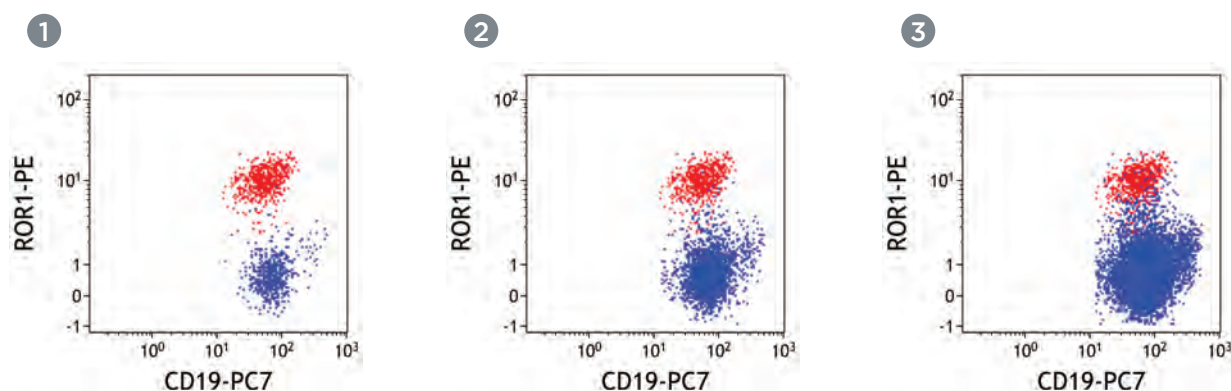

PLOT #	NORMAL B CELLS (# OF EVENTS)	ABNORMAL B CELLS (# OF EVENTS)
1	588	582
2	3,103	582
3	16,882	582

Data was acquired from a whole blood sample using a Navios Cytometer. Kaluza 1.5a analysis software was used to create digitally mixed data sets with different proportions of normal and abnormal B cells.

Antigen expression patterns of greatly abundant normal populations may overlap with the expression patterns of rare abnormal cell populations. To assure discrimination of abnormal from normal cells a carefully selected combination of antibodies is needed.

# DuraClone RE PANELS

## Antibody combinations for rare event detection in clinical research of blood disorders

The DuraClone RE Tubes are dry pre-formulated antibody panels optimized for the identification and characterization of rare abnormal blood cell populations in human whole blood or bone marrow.

- Pre-formulated 7-8 color antibody panels
- Unitized dry Beckman Coulter proprietary format
- Expert-proven marker combinations
- 25 tests/package including 3 compensation kits

Product	405 nm		488 nm					633 nm				Quality standard	
	PB	KrO	FITC	PE	ECD	PC5.5	PC7	APC	AF647	AF700	APC-AF700		APC-AF750
RE CLB Tube B80393 (25 tests RUO)	CD20	CD45	CD81	ROR-1	-	CD79b	CD19	CD5	-	-	-	CD43	ISO 9001 - 2008
RE PC Tube B80394 (25 tests RUO)	CD38	CD45	CD81	CD27	-	CD19	CD200	CD138	-	-	-	CD56	
RE ALB Tube C00163 (25 tests RUO)	-	CD45	CD58	-	CD34	CD10	CD19	-	-	-	CD38	CD20	

Usage of identical tandem dye-labeled antibodies in the compensation kits and the assay tubes assures accurate and easy compensation.



PB: Pacific Blue\*  
 KrO: Krome Orange  
 AF647: Alexa Fluor\* 647  
 AF700: Alexa Fluor\* 700  
 AF750: Alexa Fluor\* 750  
 APC-AF700: APC-Alexa Fluor\* 700  
 APC-AF750: APC-Alexa Fluor\* 750

# HIGHLY DOSED ANTIBODY COMBINATIONS TO FIND THE FEW AMONG THE MILLIONS

The achievable sensitivity in rare event analysis is mainly determined by the acquired cell number. For the DuraClone RE panels we have worked with leading expert labs to assure both, sensitivity and specificity, through state-of-the-art panel configurations with elevated antibody dosing.

DURACLONE PANEL	TARGET POPULATION	PANEL FEATURES	AMOUNT OF SAMPLE / TEST
RE CLB	Abnormal rare CD5+ mature B cells	Inclusion of ROR-1 <sup>(1)</sup> , 8 colors	300 µL of whole blood
RE PC	Abnormal rare plasma cells	Inclusion of CD200 <sup>(2)</sup> , 8 colors	Up to 5x10 <sup>6</sup> leucocytes in 100-200 µL suspension
RE ALB	Abnormal rare immature B cells	Maximal drop-in options by open PE and APC slots, 7 colors	Up to 2x10 <sup>6</sup> leucocytes in 100-200 µL suspension

1. Orphan receptor tyrosine kinases ROR1 and ROR2 in hematological malignancies. Daneshmanesh AH, Porwit A, Hojjat-Farsangi M, Jeddi-Tehrani M, Tamm KP, Grandér D, Lehmann S, Norin S, Shokri F, Rabbani H, Mellstedt H, Österborg A. *Leuk Lymphoma*. 2013 Apr;54(4):843-50.

2. Expression profiles of novel cell surface molecules on B-cell subsets and plasma cells as analyzed by flow cytometry. Llinàs L, Lázaro A, de Salort J, Matesanz-Isabel J, Sintés J, Engel P. *Immunol Lett*. 2011 Jan 30;134(2):113-21.


# STREAMLINE YOUR WORKFLOW WITH DuraClone RE

---


In the DuraClone RE kits every assay tube contains a layer with the dried-down antibody panel. The elimination of antibody pipetting shortens the turn-around-time, minimizes human error sources and purges the workflow from repetitive actions that are expensive to automate.



## Workflow For RE PC And RE ALB Panels

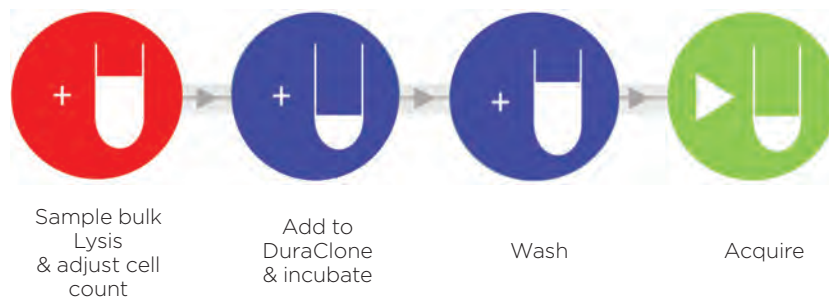


## Workflow For The RE CLB Panel

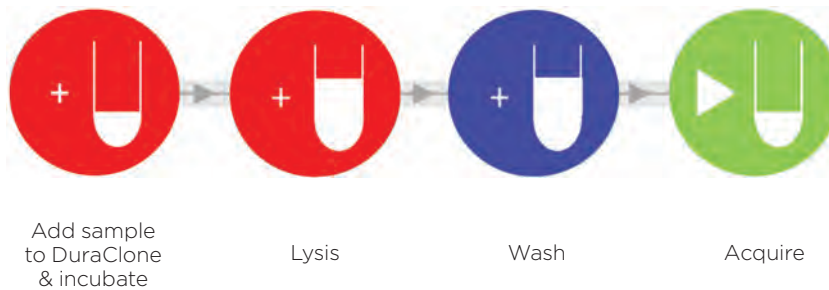



# DuraClone RE DATA EXAMPLES

## RE CLB

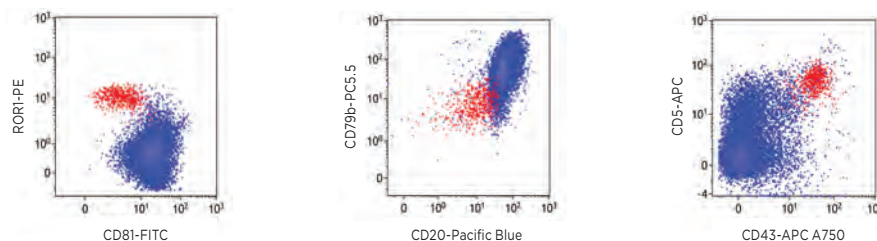


16,742 normal B cells  
577 abnormal B cells  
= **0.070%**  
of CD45+ (826,000)

### Detection of CD45<sup>+</sup>CD19<sup>+</sup>ROR-1<sup>+</sup>CD5<sup>++</sup>CD43<sup>++</sup>CD81<sup>dim</sup>CD20<sup>dim</sup>CD79b<sup>dim</sup> B cells

DuraClone RE CLB Tube	PB	KrO	FITC	PE	ECD	PC5.5	PC7	APC	APC-AF700	APC-AF750	Quality Standard
B80393 (25 tests RUO)	CD20	CD45	CD81	ROR-1	-	CD79b	CD19	CD5	-	CD43	ISO 9001-2008

## RE PC

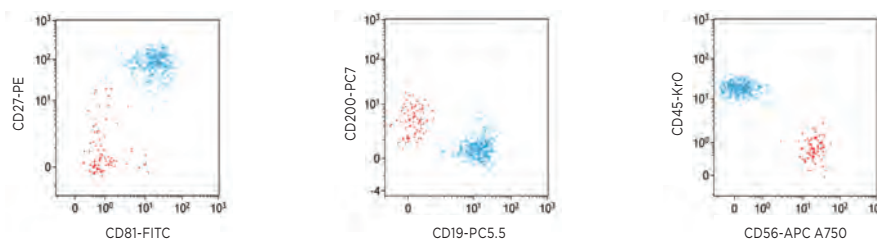


380 normal plasma cells  
84 abnormal plasma cells  
= **0.004%**  
of CD45+ (2,099,000)

### Detection of CD45<sup>dim</sup>CD38<sup>+</sup>CD138<sup>+</sup>CD56<sup>+</sup>CD200<sup>+</sup>CD19<sup>dim</sup>CD27<sup>dim</sup>CD81<sup>dim</sup> plasma cells

DuraClone RE PC Tube	PB	KrO	FITC	PE	ECD	PC5.5	PC7	APC	APC-AF700	APC-AF750	Quality Standard
B80394 (25 tests RUO)	CD38	CD45	CD81	CD27	-	CD19	CD200	CD138	-	CD56	ISO 9001-2008

## RE ALB

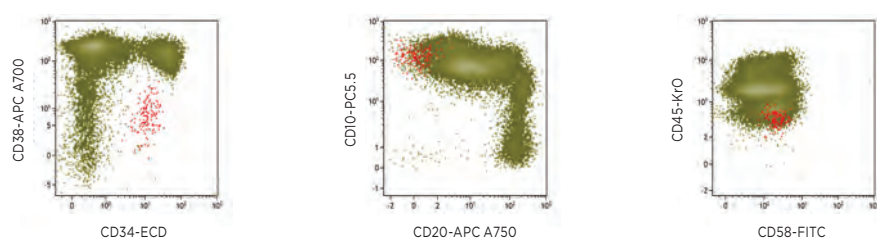

39,898 normal B cells  
132 abnormal B cells  
= **0.011%**  
of CD45+ (1,228,000)  
(merged data sets)

### Detection of immature B cells with abnormal pattern of CD10, CD20, CD34, CD38, CD58

DuraClone RE ALB Tube	PB	KrO	FITC	PE	ECD	PC5.5	PC7	APC	APC-AF700	APC-AF750	Quality Standard
C00163 (25 tests RUO)	-	CD45	CD58	-	CD34	CD10	CD19	-	CD38	CD20	ISO 9001-2008

Open PE and APC slots for maximal drop-in choice

# MORE DuraClone PANELS

## DuraClone IM for Immune System Research

The DuraClone IM antibody panels provide a comprehensive image of the immune system applying dry expert-proven antibody combinations that are easy to implement.

Product	405 nm		488 nm					633 nm				Quality Standard	
	PB	KrO	FITC	PE	ECD	PC5.5	PC7	APC	AF647	AF700	APC-AF700		APC-AF750
IM Phenotyping Basic Tube B53309 (25 tests RUO)	-	CD45	CD16	CD56	CD19	-	CD14	CD4	-	CD8	-	CD3	ISO 9001-2008
IM B Cell Tube B53318 (25 tests RUO)	IgM	CD45	IgD	CD21	CD19	-	CD27	CD24	-	-	-	CD38	
IM T Cell Subsets Tube B53328 (25 tests RUO)	CD57	CD45	CD45RA	CCR7	CD28	PD1	CD27	CD4	-	CD8	-	CD3	
IM Dendritic Cells Tube B53351 (25 tests RUO)	HLA-DR	CD45	CD16	Lineage**	-	CD1c	CD11c	Clec9A	-	-	CD123	-	
IM TCRs Tube B53340 (25 tests RUO)	TCRV $\delta$ 2	CD45	TCR $\gamma$ $\delta$	TCR $\alpha\beta$	HLA-DR	-	TCRV $\delta$ 1	CD4	-	CD8	-	CD3	
IM Treg Tube B53346 (25 tests RUO)	Helios	CD45	CD45RA	CD25	-	CD39	CD4	-	FoxP3	-	-	CD3	
IM Granulocytes Tube B88651 (25 tests RUO)	CD15	CD45	CD294	-	CD16	CD33	CD11b	PD-L1	-	-	Lineage***	CD62L	
IM Count Tube C00162 (25 tests RUO)	-	-	CD45	Counting Beads	-	7-AAD	-	-	-	-	-	-	

\*\* : CD3 / CD19 / CD20 / CD14 / CD56

\*\*\* : CD3 / CD14 / CD19 / CD56

## DuraClone IF For Immune Function

The DuraClone IF antibody panels reveal immune function at the single cell level utilizing a streamlined and simplified intracellular staining workflow.

Product	405 nm		488 nm					633 nm			Quality Standard
	PB	KrO	FITC	PE	ECD	PC5.5	PC7	APC	AF700	AF750	
IF T Activation B88649 (25 tests RUO)	CD4	-	IFN $\gamma$	TNF $\alpha$	-	-	IL-2	-	CD8	CD3	ISO 9001-2008
IF T Helper Cell (2017)	IL-17	-	IFN $\gamma$	-	-	-	IL-4	CD4	-	CD3	

## Choose Beckman Coulter for Benchmark Expertise and Innovation

For over 80 years Beckman Coulter has driven innovation. We remain committed to shaping flow cytometry technology to fit seamlessly into your lab's workflow and to provide an optimal user experience. When you choose a Beckman Coulter product you're choosing expertise, innovation, and quality assurance.



Contact your sales representative



or visit <http://www.beckman.com/coulter-flow-cytometry/duraclone-rare-event-solutions>



\* Alexa Fluor and Pacific Blue are registered trademarks of Molecular Probes, Inc.

© 2016 Beckman Coulter, Inc. All rights reserved. Beckman Coulter, the stylized logo, and the Beckman Coulter product and service marks mentioned herein are trademarks or registered trademarks of Beckman Coulter, Inc. in the United States and other countries.

For Beckman Coulter's worldwide office locations and phone numbers, please visit "Contact Us" at [beckman.com](http://beckman.com)

FLOW-2230SB01.17